
FlexiRoll™

Operator's Manual
Manuel de l'opérateur
Bedienungshandbuch
Manual del operador
Manuale d'uso
オペレーターマニュアル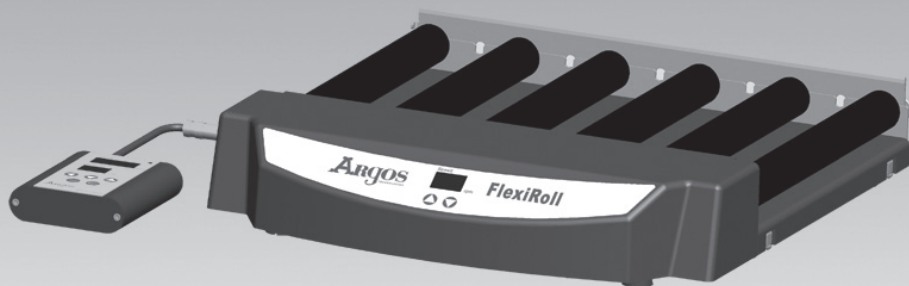

ARGOS
TECHNOLOGIES *manufacturing solutions for science*

Argos Technologies, Inc.
customer-service@argos-tech.com
Toll Free: 800-886-8675 International:
847-622-0456 www.argos-tech.com

Table of Contents

Introduction	3
Package Contents	3
Specifications	4
Important Safety Information	4
Installation	5
Installation for Stacked Units	6-7
Special Application	8
Operating Instructions	9
Maintenance and Cleaning	10
Troubleshooting	10
Warranty	11
Contact Information	11

Introduction

The Argos FlexiRoll is a digital cell roller that can be used with a variety of bottles, tubes and flasks. Applications include growth and culturing of adherent cells, suspension cells, mixing blood or vacutainer™ tubes, washing nucleic acid blots or performing hybridization protocols. The unit can be operated in temperatures ranging from 4°C - 60°C and can be used safely in atmospheric environments of up to 95% non-condensing humidity.

Package Contents

Upon receipt of the FlexiRoll, please carefully unpack your unit to ensure it is complete and in good working order. Should you find the unit damaged upon inspection, notify your freight carrier at time of receipt as well as your distributor.

Should any parts listed below be missing or you discover concealed damage once you install the FlexiRoll, contact Argos Technologies immediately for resolution (see page 11 for contact information).

FlexiRoll: 1 Unit

Cat. Number R1000 (U.S. Adapter), R1000E (Europe Adapter)

Cell Roller Rods: 6 Each

Cat. Number R1100

100-240V Power Supply: 1 Unit

Cat. Number R1101

Remote Digital Controller: 1 Unit

Cat. Number R1200

O-Rings: 2 Sets

Cat. Number R1102

Operator's Manual: 1 Unit

Warranty Card: 1 Unit

Optional Accessories:

Stacking Rods: 4 Each

Cat. Number R1205 (sold separately)

Specifications

Model	R1000	R1000E
Roller Speed	0.5 - 80 RPM	
Roller Bars Dimensions (mm)	Ø28 x 259	
Available Tubes or Bottles	Max. ø120 bottles, tubes	
Dimensions W x D x H (mm)	360 x 402 x 72	
Timing Range	0 - 999min. or continuous run	
Max. Load (Kg)	6.5	
Power Requirement	10W	
Ambient Temperature	4°C - 60°C	
Relative Humidity	<= 95%	
Net Weight (Kg)	6.5	
Power Supply	AC100V-240V, 50Hz-60Hz	
Materials	Housing: ABS Plastic / Base Plate: Stainless Steel / Rollers: Aluminum 6061	
Capacity	3 – 850cm ² Roller Bottles 20 – 15&50mL Centrifuge Tubes (requires 11 tubes)	

Important Safety Information

- Check to ensure power supply is inserted into a properly ground socket.
- Be sure to unplug the FlexiRoll when removing or installing cell rollers.
- Place the unit on a level, sturdy surface to avoid vibration.
- Never move the unit while in operation.
- Not to be used to mix flammable samples; avoid aggressive solvents or vapors.
- Do not use the FlexiRoll in hazardous, flammable or combustible environments.
- Take appropriate precautions when mixing hazardous samples.

Installation

The Argos FlexiRoll is delivered completely assembled and requires no additional set up. To accommodate the different size bottles, flasks and tubes, the cell rollers can be removed or added depending on application.

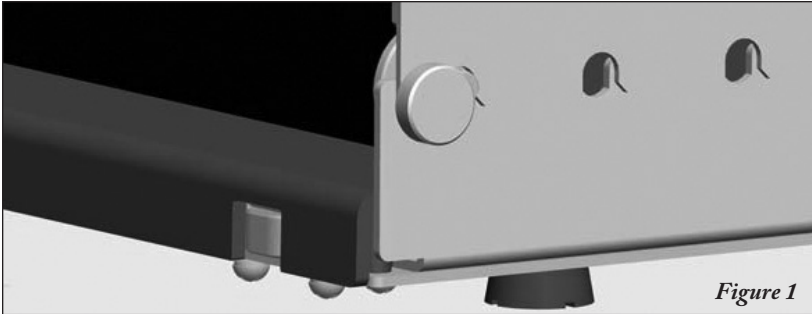
Figure 1

1. Twist the round securing knob counter-clockwise until it becomes loose. It does not need to be removed entirely. (Figure-1)


Figure 2

2. Lift up the metal retaining strip which will allow you to remove one or more of the cell rollers.
3. To install the cell rollers, follow instructions 1 & 2 above to lift the top of the retaining strip. Align the hexagon shaped hole at the end of the cell roller so that it can insert into the corresponding hexagonal shaft located on the base unit housing (Figure-2). Lay the other end of the roller in the semi circular opening on the bottom half of the retaining strip.
4. Lower the top half of the retaining strip and twist the round screwing knob clockwise to secure the retaining strip and cell rollers.


Installation for Stacked Units

Two FlexiRoll units may be stacked to maximize limited space in incubators on the laboratory bench.

Note: Stainless steel stacking rods must be purchased separately.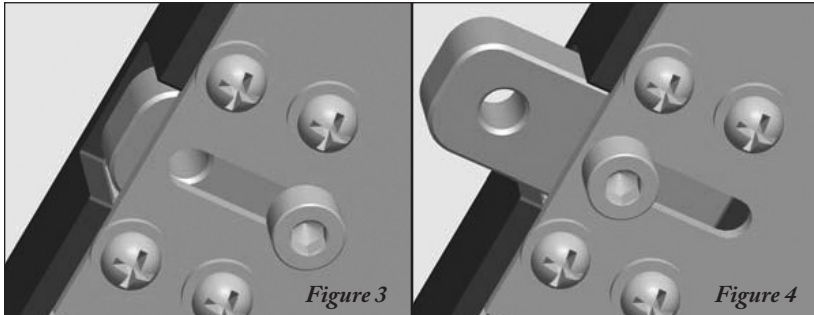

1. Unscrew the hexagonal screw on the bottom of the unit (Figure-3) until you are able to slide out the support base from the unit housing (Figure-4). Retighten the hexagonal screw to lock in place (wrench provided).
2. Repeat step #1 until all four support bases are extended and locked in position.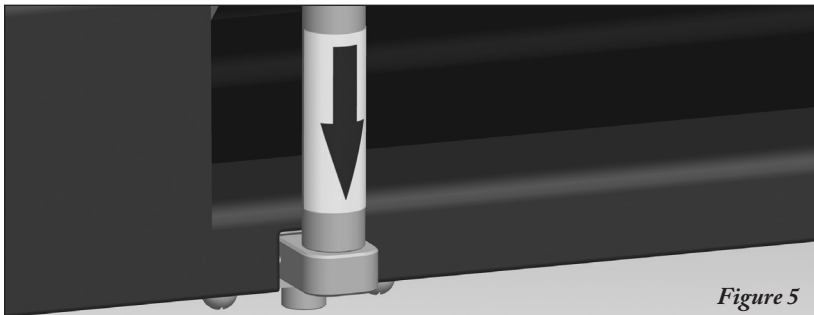

3. Insert stacking rod into the bottom support base and secure with the screw provided (Figure-5).
4. Repeat step #3 on the other stacking rods, securing in the same manner.
5. On the second FlexiRoll unit, repeat steps 1 & 2 to expose the support bases from the unit housing.

Installation for Stacked Units (cont.)

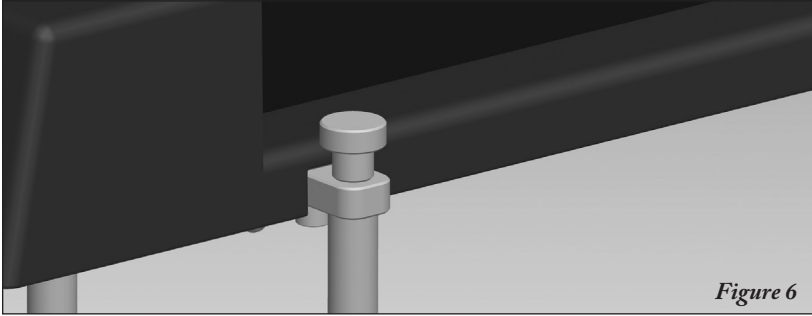
Figure 6

6. Carefully place the second unit on top of the first by setting the units support bases on top of the stacking rods. Secure with the screws provided (Figure-6).

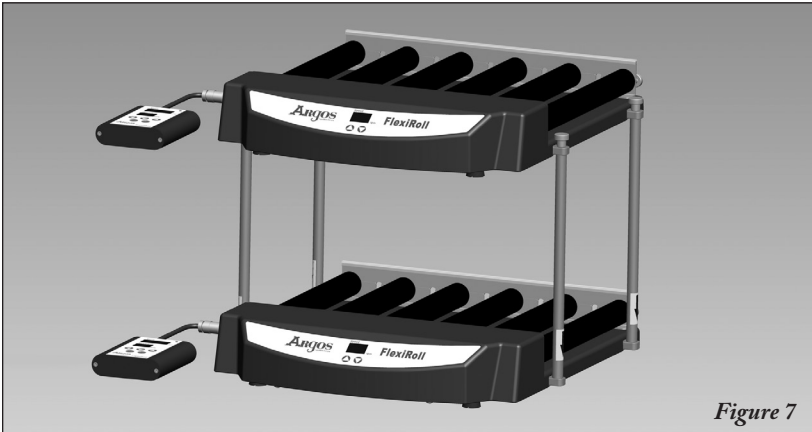
Figure 7

7. Fasten the cap to the top of the stacking rod (Figure-7).

The FlexiRoll should look like the unit in Figure-7.

Special Application

English

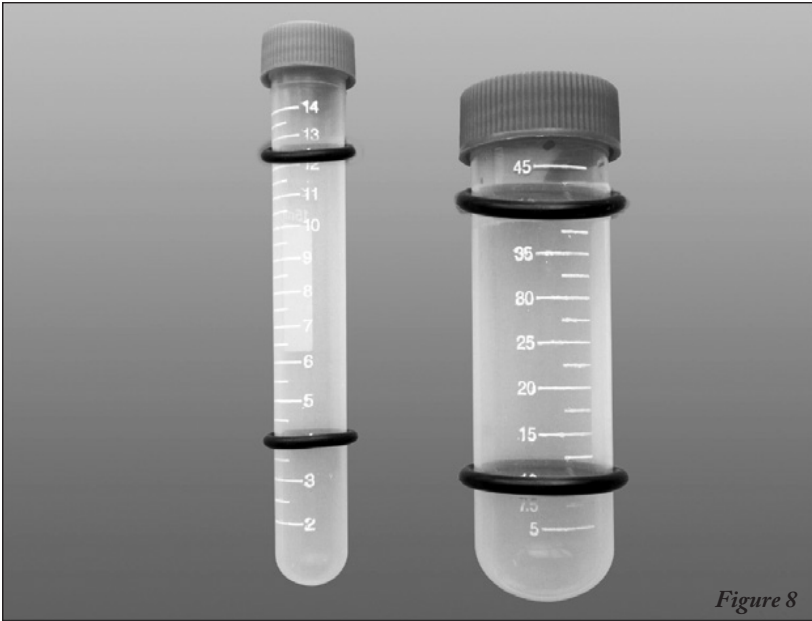
Figure 8


Special Application: 15&50mL Tubes

Please note that when rolling 15&50mL tubes, you will need to place 2- o-rings (provided) around the bottom of the tube (Figure-8). This is to keep the tubes stable in the cell roller, offsetting the larger diameter of the tube cap.

Operating Instructions

The FlexiRoll can be operated in two modes (Front Panel or Remote Digital Controller):

Front Panel Operation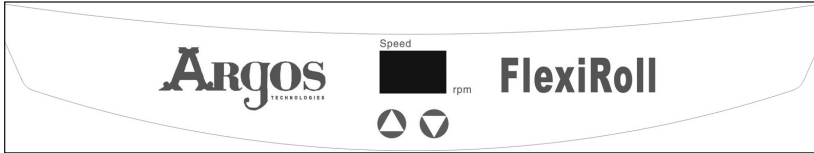

1. The FlexiRoll is activated by pressing the up arrow key to the desired speed in RPM. The lowest setting is .5 RPM and it will increase in .1 RPM increments until reaching 1 RPM. (There is no independent on / off key).
2. The speed of the FlexiRoll will be displayed in the LED panel on the front of the unit.
3. Pressing the down arrow key will reduce the speed of the unit, and is also used to turn the unit off by depressing the key to 0, the digits will flash 5 times before powering off.


Or to immediately shut off the unit, press the up and down keys simultaneously until the digital display powers off.

Remote Digital Controller Operation

In applications where the FlexiRoll is used inside an incubator, you may wish to use the remote digital controller. This allows the operator to control the time and speed of the FlexiRoll.

Note: When utilizing the remote digital controller, the LED display on the unit will not operate.

1. Insert remote digital controller cable to the socket located next to the power supply.
2. Depress "set" key once to activate time setting. The flashing digit can be adjusted by depressing the up or down key. To move the cursor to next digit, press the left arrow. This will allow you to set the time up to 999 minutes.
3. To set the speed, once again depress the set key and the speed digits will flash. Using the up, down or side arrow key will allow you to set the FlexiRoll to the desired speed.
4. Press the start key once to start operation. By pressing a second time, the unit will stop.
5. To rapidly increase / decrease the speed and or time settings, depress the up or down arrow key for 3 seconds.

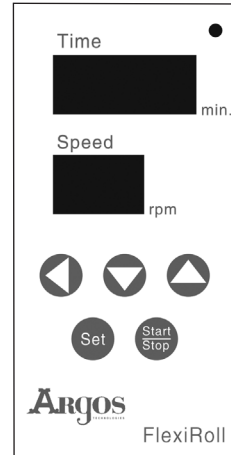Remote Digital Controller

Note: The timing function is deactivated when the unit is used in continuous mode (reads "00" in time display).

Maintenance & Cleaning

- No lubrication is required.
- Before cleaning be sure to unplug the power cord from the wall socket.
- The housing can be cleaned using a soft damp cloth and mild non-corrosive detergent.
- After cleaning, ensure that all parts are thoroughly dry before attempting to operate the unit.
- Do not immerse the FlexiRoll in liquid.

Troubleshooting

Error	Cause	Solution
Unit does not start	Power cable not connected.	Check that power cable is firmly plugged into wall outlet.
	Power Supply.	Check that there is power to the wall outlet.
	Remote Control.	Check that cable is correctly installed into the unit.
	Sample Overload.	Ensure total weight does not exceed 6.5Kg.
No Display	Power Failure.	Ensure unit is powered on.
	Timer cycle completed.	Repeat run.
LED Light On; unit not working	Overloaded.	Check and repeat run.
Non-responsive Keypad	Remote digital controller has been connected.	Detach remote digital controller from unit.

Warranty

Argos FlexiRoll is warranted to be free of defect in materials and workmanship for a period of 12 months from the date of purchase, providing that:

- The unit has not been opened or attempted to be opened by the end user.
- The user ensures that it has been carefully cleaned and decontaminated.
- The unit has been operated according to the instructions and not in any improper way.

In order to qualify for warranty repair; the customer must return the warranty card supplied with the unit that includes all of the information required to validate proof of purchase. Should warranty service be required, repair or replacement will be provided by Argos without charge. Freight shipping the unit to Argos will be the responsibility of the purchaser. There are no other warranties or guarantees, implied or otherwise expressed including warranties of application or merchantability.

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